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1	Main Design Data	<p>ReactionFlask (L) : 100, FlaskNicks: 5 or 6, Stirring Power: 370W, RotationSpeed: 50–900rpm, Torque: 22000g. cm, ElectricalRequirements: 220V, 50Hz.</p> <p><b>Material: 304 Stainless frame.</b>  <b>Standard configuration:</b></p> <ol style="list-style-type: none"> <li>1. Capacity – 100L      Answer: yes</li> <li>2. Operate temperature – 70 °C      Answer: no problem</li> <li>3. Operate pressure – atmospheric      Answer: Yes</li> <li>4. Rotation speed – 500 rpm      Answer: no problem.</li> <li>5. Material of stirrer’s shaft: - 304 with PTFE cover      Answer: No problem</li> <li>6. Motor: explosion-proof      Answer: No problem</li> <li>7. Frame material: 304      Answer: No problem</li> <li>8. Flask Necks – 5      Answer: No problem.</li> </ol> <ol style="list-style-type: none"> <li>1. Constant pressure funnel: 5 l</li> <li>2. Condenser with bottom receiving flask</li> <li>3. Thermocouple</li> <li>4. big feeding opening (D = 50 mm) with PTFE cover</li> <li>5. Vacuum Mouth</li> </ol> <ol style="list-style-type: none"> <li>9. Could you specify the size for each neck? Pay attention, please, the condenser no need! But we ask the addition funnel instead of it.</li> </ol> <p>Answer: for the size of each neck: now we will tell following:  Middle: 60mm, Left is 40mm, Right 40mm, Behind(rear): 34mm,</p>
---	------------------	--

		<p>Ahead:50mm.</p> <p>10. What size of nozzle for connecting water tube to jacket? Could you specify it in your quotation?</p> <p>Answer: outer Diameter: 8mm.</p> <p>11. Metallic protection shield is not needed      Answer: Ok.</p> <p>12. Distance from floor to the drain valve shall be 60 mm Answer: ok.</p> <p>13. In addition to the reactor, the supply should include:</p> <ul style="list-style-type: none"> <li>a. glass stopper for every neck (3 pieces for each neck )</li> <li>b. The PTFE cover for big opening (3 pieces)</li> <li>c. O-ring (F-rubber) for flush bottom valve, 10 pieces. Necessarily check them for resistance to acids</li> <li>d. Spare flush bottom valve (1 pieces)</li> <li>e. Constant pressure funnel - (4 pieces)</li> <li>f. Vacuum Mouth – (4 pieces)</li> </ul>
2	Discription	<p>Circulation liquid of hot, cold or cryogenic can be filled into the jacket of reactor. Substance reacts by stirring under normal or negative pressure in aieproof glass container. Reaction process can be watched clearly through crystal reaction flask. Jacket Glass Reactor is ideal equipment in library experiment or medium-production in moden biology and pharmcy, and is also applied in synthesizing materials in fine chemical industry. Features in Design: 1. Integrated jacket reaction flask has big caliber with good seal and high strength. 2. Professional PTFE cover supports high speed stirring in vacuum. 3. Three standard mouths can be used for umlti-purpose according to customer needs. Standard fitting are chatging cork, thermometer tube, and reflux condenser. (Customization available for special needs) 4. More designs on safe and easy operation are applied on models of 20L and above.</p>
3	Note	<p>max speed: 500rpm</p>

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Thank you for using HEB brand jacketed glass reactor, our motor is good service and good quality. If you have any technical questions, plz contact with our technician, Tel:86-29-81047996

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Mainly describe how install and use for jacketed glass reactor.

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The misunderstanding of operation will result some accident, reduce product life and function of product, so plz read our instruction before use.

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To give final customer for our instruction.

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Plz put instruction in good postion so that read it any time.

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Only is reference for our instruction.

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Must pay attention to item following, for good operation.

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Plz must read all contents of instruction before  
intall, use, maintain, examine.

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The electrical requirement must be conformance with our  
instrument precribed.

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To prohibit high speed rotation when it is no load.

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Plz keep speed controller and motor drying

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Cut off power inside 5mins,plz pay attention to security by  
discharge of speed controller and capacity.

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Plz pay attention to hard thing(watch, ring, etc) to prevent from damage of crash

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Must be careful for connecting between water pipe and vacuum valve or deflation valve(rear of lid) to prevent glass crack result in people wound.

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Plz attention to security when use heating circulation system, we suggest use heating oil of small dangerousness(for instance two-A silicon oil).plz pre-open circulating pump when user is going to get circulation, slowly rising temperature under circulation of heating oil.

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### Summery

The jacketed glass reactor often be used as bio-chemical instrument, it is used in refined chemicals, pharmaceutical, scientific experient, etc, user can do concentration, distillation, seperation, extraction under constant speed, constant temperature.

For your convenience, jacketed glass reactor, the fully assembled reactor head, stirrer shaft or agitator, stirrer shaft coupling, Tflon stirrer bearing, temperature contoller and speed controller, s. s inlet/outlet adaptors, and motor are shipped assembled on support frame.

Unpack all parts and check against packing slip to make sure you have received all necessary components, if it is possible, plz keep some of packing material so that you can use it continously, to reture us to repair our machine.

Cratel: Reactor Vessel(plz be carefully when move it), assembled motor, assembled stirrer shaft or agitator,

500ml receiving flask with three joints, all screws, all  
“O” ring and half “O” ring, Glass Lid, Condenser, Vacuum  
Valve, wrenches, charging valve, discharging valve, etc.

s. s. frame. plz install them carefully to reach good balance.

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Sealing for using PTFE materials and special rubber combined  
sealing ring to endure various kind of different solvents and  
keep high vacuum.

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High efficiency condenser to make sure high receiving rate.

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Standard connection way to unload easily.

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To install discharging switch, it is quite easy for receiving  
hadralytic enzymes.

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To install vacuum meter, to choose the best vacuum under  
materials of low boiled point.

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The structure is reasonable, main parts were made of high quality stainless, aluminium alloy and other new sythintic material, its priority includes resisting corrosion, dust, and enduring various kind of solvent.

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Gls parts were made of GG17 silicon glass material, to own good physic and chemical function.

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Electrical parts were made of high quality processed part imported.

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Adjusting speed is no class for speed controller, motor no flame, it is reliable and secure, long life(to choose imported precessed speed controller above 30L).

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After open package, check up all parts

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Install stainless frame to use stainless pipe and metal fixing.

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Instll speed controller on right rod of rear(plz look at picture), then insert plug; vacuum meter will be installed left of rear, then firm it well.

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To firm “0” ring on frame(plz look at picture)according to your desired height, and put flask vessel on “0” ring;to firm neck of flask vessel to use half” 0” ring, total two half” 0” ring, to screw up after fold two half” 0” ring, plz pay attention to verticality of flask vessel when you install.

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Stirrer is firmed on head of “gimbal joint”, then stirrer inserts into hole of Column from center of glass lid

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The slipper of rear rod is to adjust different direction, it is a kind of clamp.

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The right side of lid for 40#standard joint with condenser, the lift side of lid for 40#standard joint with charging funnel, ahead of lid for 24#standard joint with pipe of Phmeter, rear of lid for 34standard joint, it is a multi-function joint.

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Must keep cleaning when user install glass parts, because of every joint with grease oil.

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To press gimbal joint to make stirring, if its stiring is steady, the adjustment is ok for this stirring rod.

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Check up electrical requirement is conformance with machine required.

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To power on, open switch of power from speed controller, then choose desired speed to use speed botton.

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Flowing of material and rotating speed of motor could occur(sympathetic vibration), then plz change speed of motor to avoid it.

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Must rotate to use hand after stirrer is installed, look at if it is vertical, if it is not vertical, plz adjust it to reach good verticality, then open power, increasingly adjust speed from slow to fast.

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if you need to adjust stirrer, plz adjust up or down about 5cm.

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If user find sealing is not good, plz check up rotary axletree of jacket of stirrer.

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All parts are made of GG-17 material, plz make sure glass material is good.

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Outside of jacket with circulation joint, user can do heating reaction to infuse heating oil, can do cooling reaction to infuse cooling liquid, if user want to use heating of 蒸汽, it is not overcome 0.5K-1K. (note: we suggest to do not heating of 蒸汽)。

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Discharging valve of bottome will get fox when user do low temperature, plz unfreeze when user uses discharging valve.

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If user find many granules in liquid, they could be acumulated in stopcock of discharging valve. plz must clean it when you are using reactor, again.

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To use soft clothing to erase all joints, then coat a little of vacuum grease.

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Don' t tightly screw on all joint, plz periodically move to avoid tight connection.

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First open switch of power, then operation from slow to fast, plz keep machine stopping after use, close switch.

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The switch of PTFE, plz don' t strongly screw on it in case glass is damaged

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Plz eresease various kinds of oil. durty, solvent residual after use.

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Periodically, plz clean some sealing ring. Method: take down sealing ring, check up if it has durty on axis, then clean it.

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Electrical part, plz prohibit humidity.

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Must purchase standard part from our factory, plz don' t randomly buy other sparepart from other factory.

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Plz cut off power or water when we check and repair machine.

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Plz maintain it after use three year(only suggestion)

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Note: our reactor is fragile product, plz carefully use it, and we don' t repair fragile product(glass), so if some customer need fragile product, need to pay cost.

## 六、故障简易排除 Malfunction Removing.

Phenomenion	Reason	Removing
开启电源开关, 指示灯不亮。The Indicator is not light when you power on	External source no pass or other	请专业电工检查电源, 插座。Plz check up power supply and jack
	Short circuit	将电源开关置于 OFF 位置, 再换置保险管。Plz put switch to OFF, exchange new fuse.
电源指示灯亮, 但不旋转。Indicator led is light, but not rotary	旋转轴生锈 rotary axis is rust	停止使用, 与供应商联系 Stop to use it, plz contact with supplier.
	电机, 电气箱故障 Motor, electrotechnical box in malfunction.	
	未连接七芯插头 no connection with seven-connector electrical male plug	重新连接七芯插头 Reconnect this plug.
Vacuum is dispealed	Check up glass is ok? Glass switch is ok?	Plz replace sealing ring and switch
有真空, 但抽不上 Has vacuum, but couldn' t reach to max vacuum	密封圈磨损, 连接真空开关泄漏 Sealing or vacuum switch plz check.	Plz replace new parts
真空时有时无 sometime is in vacuum, sometime is not in vacuum	Axis has durty, loose connection	请清除污垢检查真空表, 真空泵 plz erease durty, check up vacuum meter and vacuum pump
	真空软管老化 sofy pipe aging	请更换真空软管 plz replace it.

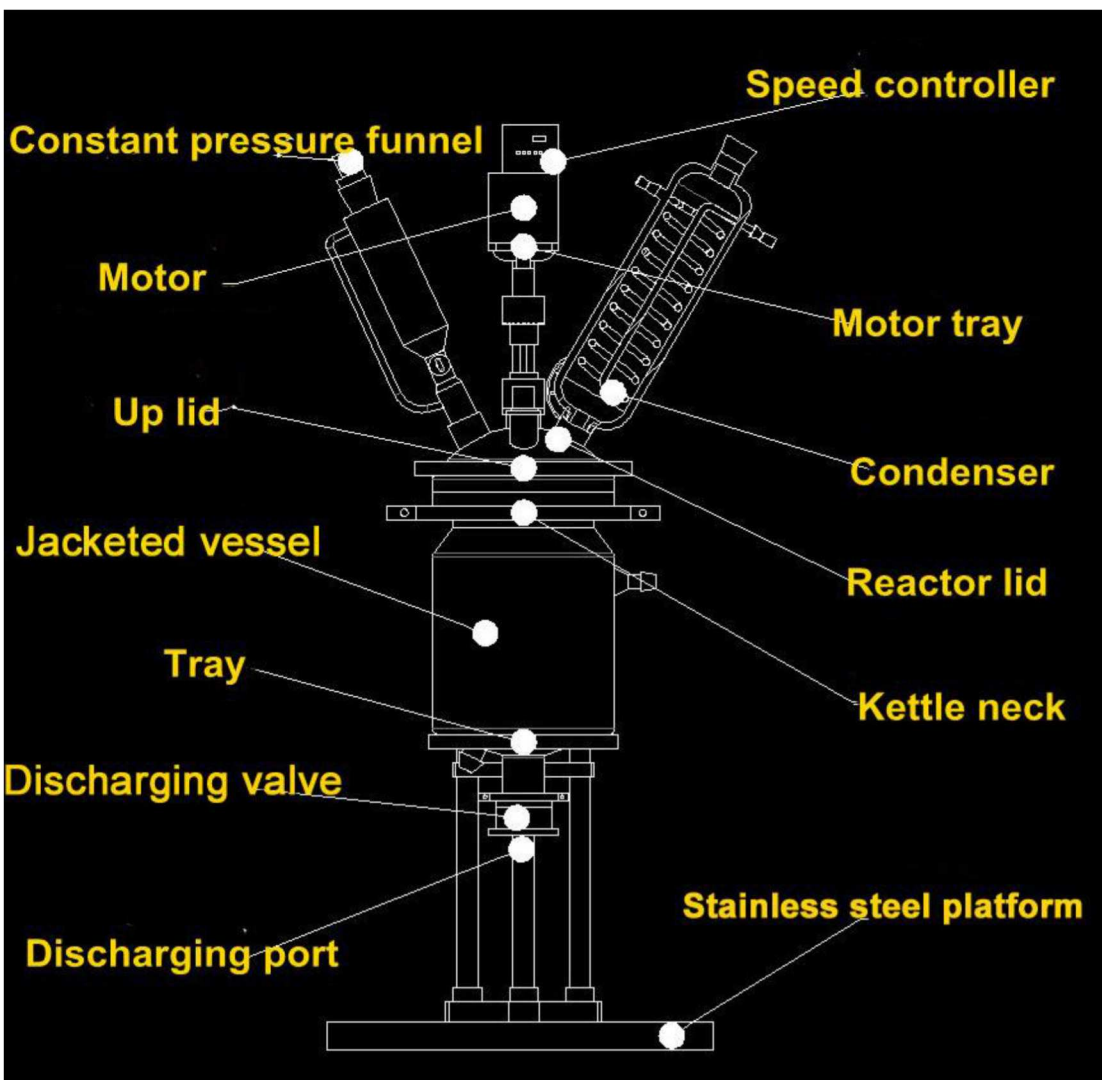
主机芯轴左右松动 Main machine	紧顶螺丝松动	旋紧
电机温度过高（室温加40℃属正常） The temperature of motor is high(RT-90degree is normal)	Overload	停机，用手使机轴转，是否很重，清除密封圈与玻璃轴接触部上的污垢，涂上真空脂 Stop working, clean up dirt between sealing ring and PTFE stirring axis
转速显示与实际不符 Speed Display is not accord with actual speed	电压不稳定，自身有误差 The voltage is not stead, or it has some error(it is normal)	与供应商联系 plz talk with supplier

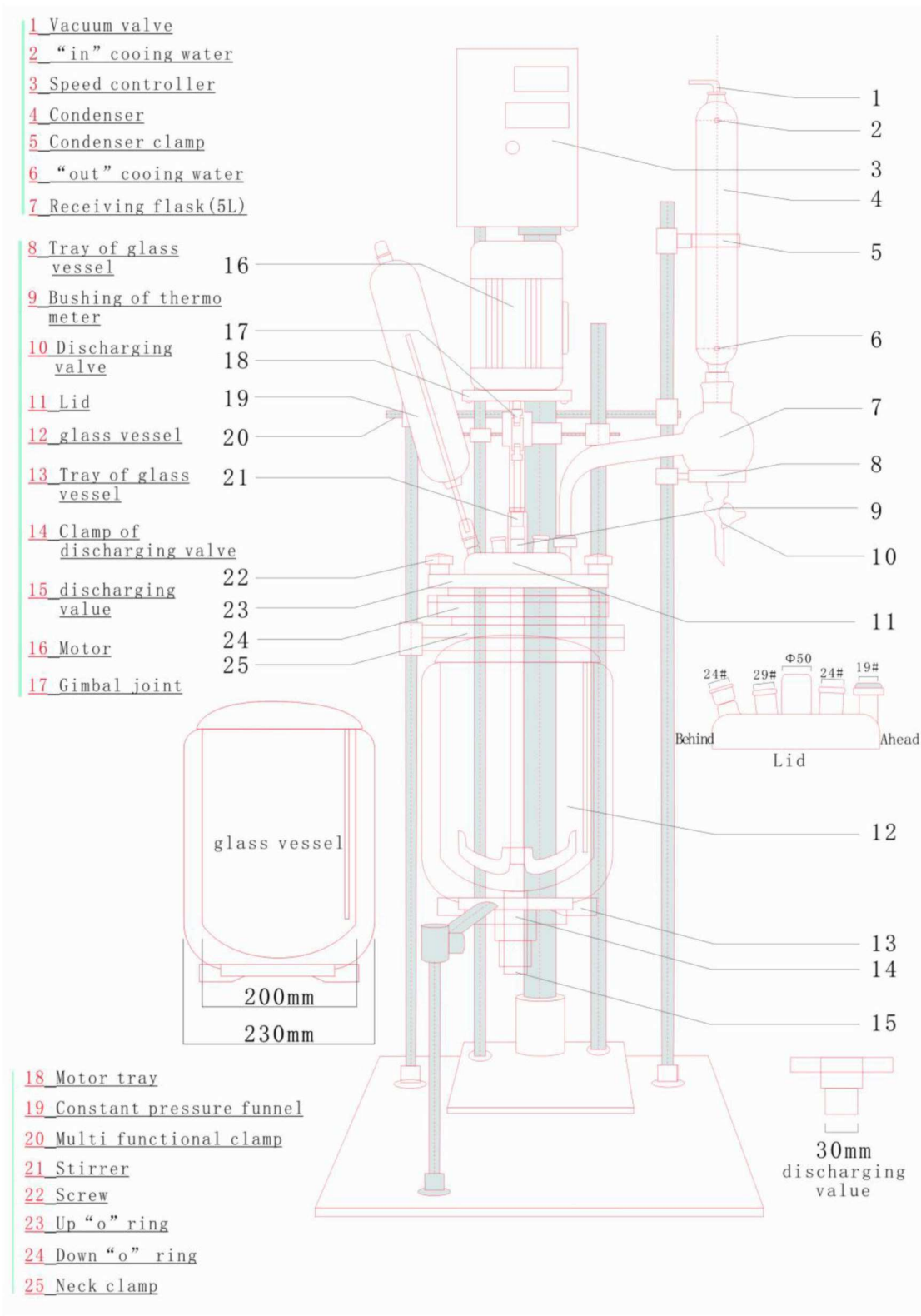
## Product Drawing(1L-5L jacketed glass reactor)

Our Part following:

- (1): HEB-FC-307: Frequency-convertor for digital display
- (2): HEB-10: Condenser
- (3): HEB-11: Electric Stirrer motor power: 40W-120W.
- (4): HEB-21: Lock Clamp
- (5): HEB: Constant Pressure Funnel
- (6): HEB-1: half-ring Clamp
- (7) HEB-07-12: receiving flask
- (8) HEB-021: black lock clamp
- (9) HEB-2: discharging Valve
- (10) HEB-99-16: 5 Neck Lid
- (11) HEB-0 Screw
- (12) HEB-RH-208 Lock Clamp "O"
- (13) HEB-5: support rod 4
- (14) HEB-RH-210 Support Ring "O"
- (15) HEB-JR-111 1L-5L Jacketed vessel
- (16) HEB-ST-101 Stirrer Shaft Stainless Teflon agitator
- (17) HEB-RH Lock 4
- (18) HEB-DS-158 Discharging Valve not existing materials
- (19) HEB-4 Main support rod
- (20) HEB-3 Flat Base

Note:


Product' s Drawing(10L-100L jacketed glass reactor)

1. Item	Part	
NO	Numbr	Description
1.	HEB-10	Condenser


2.	HEB-15	Connection Joint
3.	HEB-11	Electric Stirrer Motor(Proof-Explosive) 50-900rpm Power:370W
4.	HEB13	Pressure Meter Max pressure 0.098Mpa
5.	HEB-07-12	Receiving Flask(2L-5L Optional)
6.	HEB-99-16	5-Neck Lid diameter for Middle:60,Left:40, Right:50,Back:34mm, Front:29mm
7.	HEB-FC-307	Frequency-Convertor for Digital Display
8.	HEB-BP-502	Beaded Pipe Coupling(outlet)
9.	HEB-BP-503	Beaded Pipe Coupling(Inlet)
10.	HEB-JR-111	10-100L Jacketed Vessel
11.	HEB-ST-101	Stirrer Shaft Stainless, Teflon Agitator
12.	HEB-RH-210	Aluminium Support Ring
13.	HEB-DS-158	Discharging Valve Not Existing Materials
14.	HEB-LC-199	Heavy Duty Lockable Casters
15.	HEB-VM-201	Vacuum Mouth
16.	HEB-SF-08	Support Frame,Stainless
17.	HEB-CP-0	Constant Pressure Funnel
18.	HEB-PTFE	PTFE cover for big mouth
19.	HEB-GV	Glass Valve
20.	HEB-TP	Bushing of temperature meter
21.	HEB-FS	various kind of clamps

# JACKED GLASS REACTOR

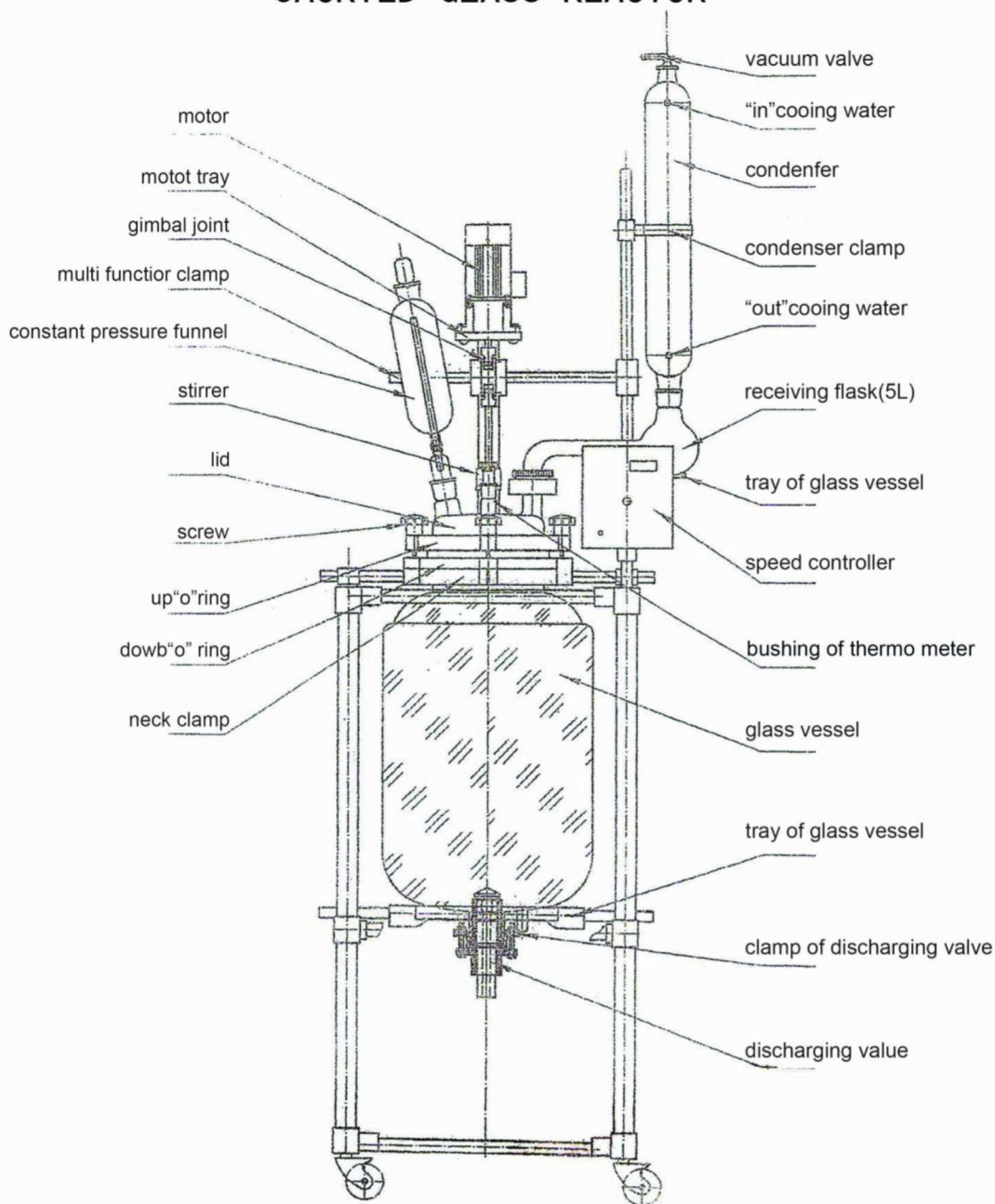Model	Parameters Inside's/outside's dia of glassvesse	Glass lid with 5 necks				
		Middle	Left	Right	Behind	Ahead
TFD-50	355*400	Φ60	40#	Φ50	34#	29#
TFD-30	310*355	Φ50	40#	40#	34#	24#
TFD-20	310*355	Φ50	40#	40#	34#	19#
TFD-10	245*260	Φ50	40#	40#	34#	19#

# JACKED GLASS REACTOR

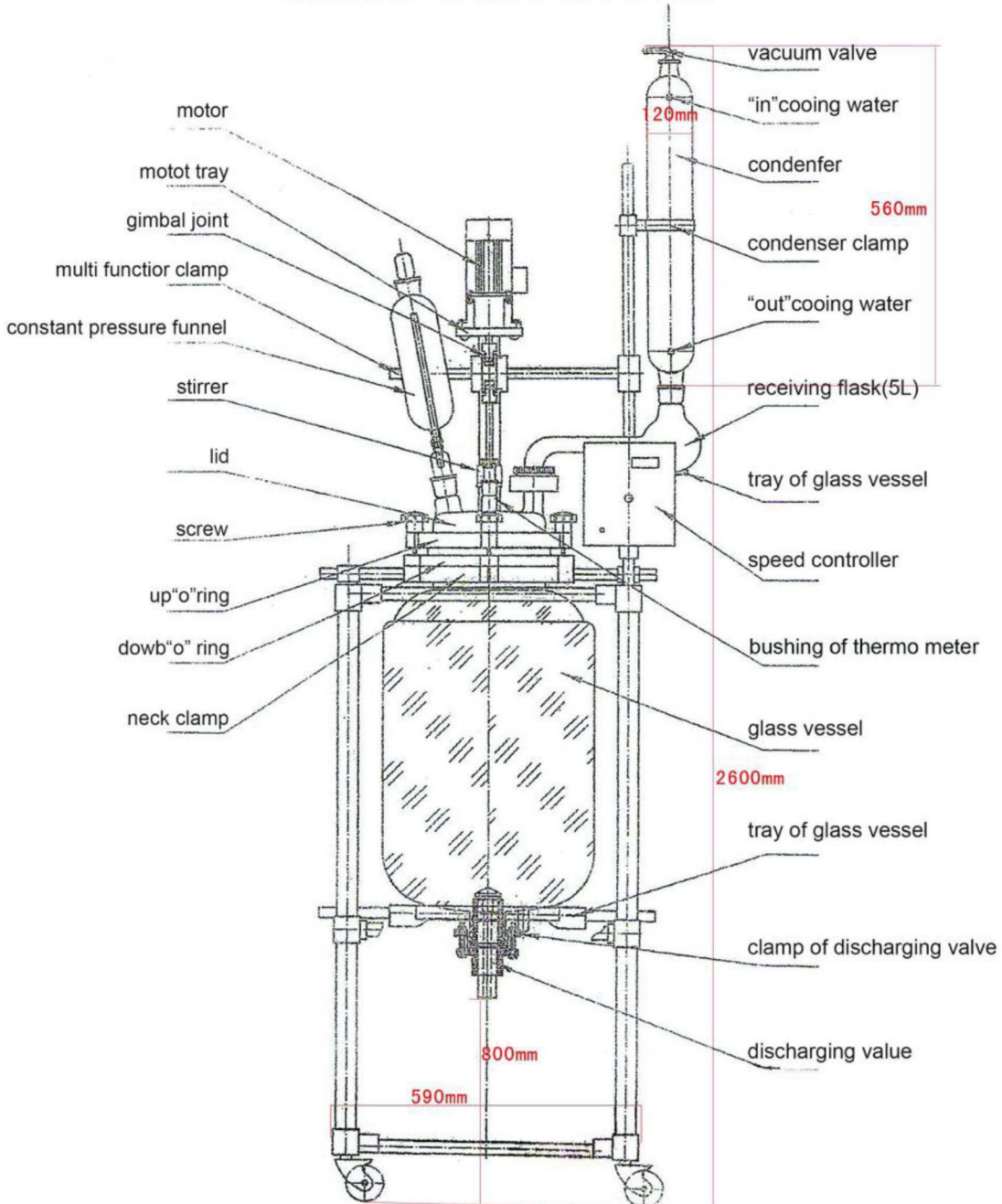Parameters Model	Inside's/outside's dia of glassvesse	Glass lid with 5 necks				
		Middle	Left	Right	Behind	Ahead
HEB-100	405*460	Φ60	40#	Φ50	34#	29#

Quality Certificate

Product Model:

Batch No:

Manufacturing Date:

Signature:

Warranty Card	<p>The product rely on Q/320211 DFX01,02-2000 standard to approve selling.</p> <p>The reason is in manufacturing, or the motor is in malfunction, warranty time for one year.</p> <p>Buying date: Invoice No:</p>	<p>Instruction:</p> <ol style="list-style-type: none"> <li>1. the user should fill in "WarrantyCard" and WarrantyRegistration Card</li> <li>2. the warranty registration card will be sent to us in one month.</li> </ol>
---------------	---	---

**Warranty Registration Card**

Company or unit			Address		
customer		Tel		P.C	
dealer		Invoice No		Buying Date	
Product model		Batch No		Mark	
Quality Tracking	Date	Repair Record			Contact person
User suggestion					

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